


Connecting a Single NIRSport2 8X8 Device


- Log into the NIRSport2 laptop and plug router into the outlet.
 - Click on *Network and Internet settings*.
 - Click on “Your access point name” broadcast.
 - If it prompts for a password, please use the following Username and Password to connect.
- Turn on the NIRSport2 machine.
 - Hold down **Power** till it turns green.




- Please confirm that the appropriate cap is populated.
- Check that all stabilizers in the cap are intact, and please replace any that are broken or loose.
- Plug in the sources, detectors, and triggers into the NIRSport2 machine.
**Note: You should hear a click when it's plugged in correctly, and please continue to keep the plugs level to the NIRSport2 machine.*

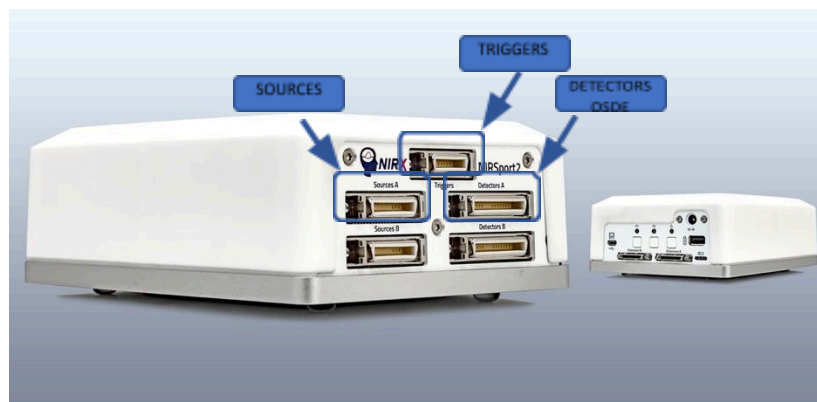


- Open **Aurora** on the fNIRS laptop.
 - Click on the 5th icon from the left-hand side of the screen. If you run your computer mouse over the icon, it should say **Open Lists of Devices**.

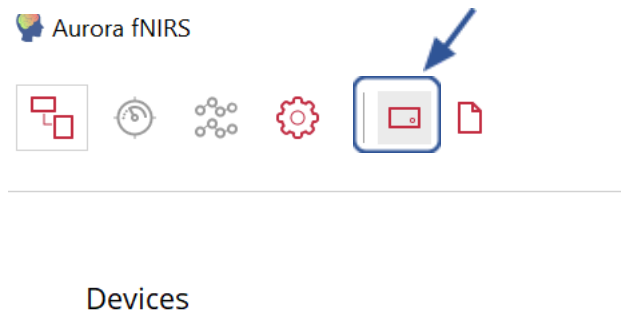

- The NIRSport2 machine ID and icon should pop up
- Then, click on **NSP2_2114_0285_A**. (this ID will change if you have purchased your own NIRSport2 device)
- If the NIRSport2 machine is on, and it does not automatically connect to the computer, please type in the Password that is located on the bottom of the NIRSport2 machine. It will prompt for it if it doesn't connect.

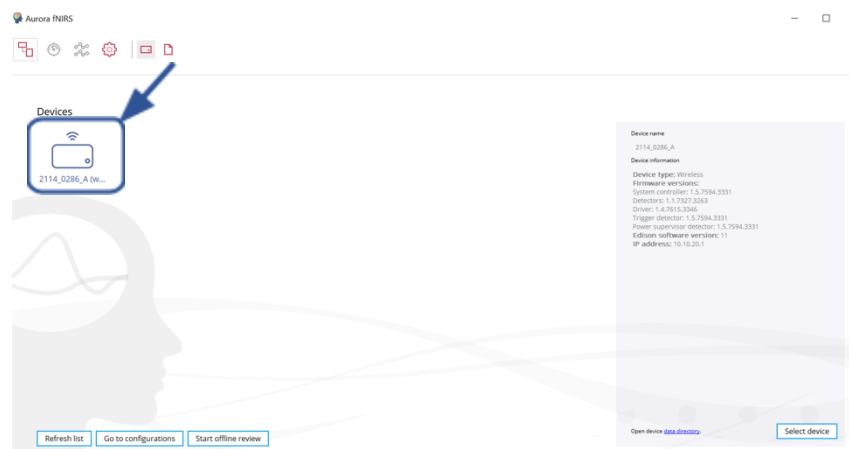

- **Note: If it does not pop up, click **Refresh List**.*

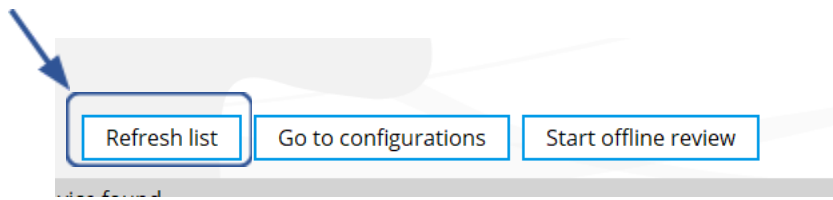

- Double-click on the NIRSport2 machine icon.
- Various configurations will pop up, please click on the configuration name for your study.
- Once selected, it will show a red checkmark and take you to a new window.