

Guide to set up FCSExpress site license

Please follow the procedures in sequence.

1. Register for De Novo Software account. Your university email is required.
<https://denovosoftware.com/register/>
2. Email XiaoxuanFan@som.umaryland.edu, include your lab name, whether you are our facility user and how many licenses your lab already has.
3. Register for site license:
<https://www.denovosoftware.com/RegisterForSiteLicense/?SignupCode=UMGCCC+FCSS>
4. You will receive the approval email within one day.
5. Download at: <https://denovosoftware.com/full-access/download-landing/>
6. Launch the installer. Choose “Stand Alone” when prompted.
7. Run FCS Express 7. Follow the wizard:
 - 1st option “Automatically through this wizard, I already have a De Novo Software account”.
 - Enter your De Novo Software Account credentials.
 - Choose configuration file for your Flow Cytometry site license.
 - Complete the wizard and FCS Express 7 will launch.