

Annexure-II

Calibration Report of Disintegration Tester (MODEL: ED2-SAPO)

Company Logo Here XX PHARMACEUTICALS LIMITED 117 Adams Street, Brooklyn, NY 11201, USA
--

Page 1 of 2

Calibration Report of Disintegration Tester (MODEL: ED2-SAPO)

Make: Model No.: Serial No.: Instrument Code No.:	Frequency: Location: Date of Calibration:
--	--

1.0 Observations:

1.1 Calibration of Temperature in Jar: (Ref. ID No. of Thermometer:)

Jar	Set / Display Temperature (°C)	Actual Temperature (°C)	Acceptance Criteria	Remarks
A	37.7° C		Display Temp. $\pm 2^{\circ}\text{C}$	
B	37.7° C			

1.2 Calibration of Timer: (Ref. ID No. of Digital Timer:)

Basket	Set Time (min.)	Actual Time (min.)	Acceptance Criteria	Remarks
A	15 min.		14.7 – 15.3 min.	
B	15 min.			

Page 2 of 2

1.3 Calibration of Stroke/min.: (Ref. ID No. of Digital Timer:)

Basket	Nominal Value (Stroke/min.)	Actual Value (Stroke/min.)	Acceptance Criteria	Remarks
A	30		29 to 32 Stroke/min.	
B	30			

1.4 Calibration of Traveling Distance: (Ref. ID No. of Scale:)

Basket	Nominal Value (mm)	Actual Value (mm)	Acceptance Criteria	Remarks
A	55		55 ± 2 mm	
B	55			

1.5 Calibration of Basket to Beaker Bottom Height: (Ref. ID No. of Scale:)

Basket	Nominal Value (mm)	Actual Value (mm)	Acceptance Criteria	Remarks
A	25		25 ± 2 mm	
B	25			

Conclusion:

Next calibration due on:

Calibration done by:

Checked by:

Approved by: