


Tata cara pembuatan SSN name (untuk Sample Set)

Jika saat akan pembuatan sample set, ternyata sample set yang akan dijalankan tidak tersedia? Missal tidak ada UV,PDA,Arc dll silahkan pake cara ini yah

- 1) Buka project empower lokasi yang akan di run
- 2) Buka pada bagian “methods”

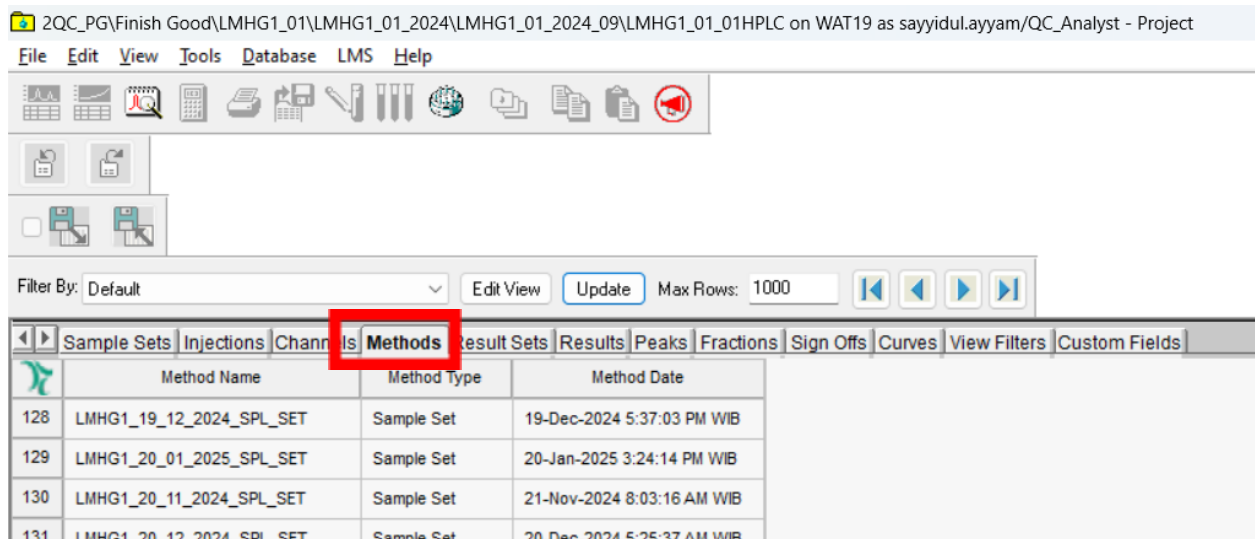

- 3) Kemudian cari contoh ssm yang sudah ada (ini contoh pada project LMHG1 hanya ada yg PDA tidak ada UV)

143	LMHG1_30_11_2024_SPL_SET	Sample Set	02-Dec-2024 8:08:57 PM WIB
144	LMHG1_31_07_2025_SPL_SET	Sample Set	31-Jul-2025 11:19:02 AM WIB
145	LMS_LMHG1_PDA	Sample Set	13-Jan-2026 3:12:00 PM WIB
146	PERSIAPAN KOLOM	Sample Set	05-Aug-2025 8:55:04 AM WIB
147	PERSIAPAN KOLOM PDA	Sample Set	30-Sep-2025 8:54:33 AM WIB

4) Open sample set contoh tersebut

Function	Method Set / Report or Export Method	R
Condition Column	MA_LMHG1_01_M01_PDA	
Inject Standards	MA_LMHG1_01_M01_PDA	
Inject Samples	MA_LMHG1_01_M01_PDA	
Inject Samples	MA_LMHG1_01_M01_PDA	
Inject Samples	MA_LMHG1_01_M01_PDA	
Condition Column	MA_LMHG1_01_M01_PDA	
Condition Column	H2O_Mtd_Set_PDA	

5) Silahkan ganti method set nya sesuai yang di butuhkan (missal diganti jadi UV pada kasus ini)

Function	Method Set / Report or Export Method
Condition Column	Mixagrip HG_LMHG1 01_01HPLC UV
Inject Standards	Mixagrip HG_LMHG1 01_01HPLC UV
Inject Samples	Mixagrip HG_LMHG1 01_01HPLC UV
Inject Samples	Mixagrip HG_LMHG1 01_01HPLC UV
Inject Samples	Mixagrip HG_LMHG1 01_01HPLC UV
Condition Column	Mixagrip HG_LMHG1 01_01HPLC UV
Condition Column	H2O_MTD_SET

6) Kemudian save as dengan nama baru yah

