

How to order DNA sequence

20161206 Uda

Use Fasmac

<https://www.bio.fasmac.co.jp/FasmacWebSystem/ja-JP/Account/Login.mvc>

Mail address: k-aoki@nibb.ac.jp

Password: okazaki

Choose 「DNAシーケンス解析」 and fill in sample name

bp of DNA is not essential (about 700bp?).

Input your e-mail address to 「追加メールアドレス」.

After input and order, **please print “order sheet” to enclose.**

Sample Preparation

Plasmid 300-600ng*

Primer (50uM) 0.13uL*2

DDW Up to 14uL

* 150ng plasmid is enough to analyze?

*2 When you use M13 primer (0.1ug/uL $\approx$ 20uM), please add 0.32uL

Write sample number to PCR tube like below



Assemble sample box and put sample tube and **order sheet** in it.

Paste delivery order sheet and take it to office at 1F.

(You can receive analysis data next day when you submit until 3p.m.).

After receiving analysis data, print e-mail [解析完了のご報告<ファスマック>] and submit it to Onoda-san.

If you have a question, please read

<https://www.bio.fasmac.co.jp/FasmacWebSystem/Download/%E3%82%B5%E3%83%B3%E3%83%97%E3%83%AB%E8%AA%BF%E8%A3%BD%E3%81%AE%E3%82%AC%E3%82%A4%E3%83%89%E3%83%A9%E3%82%A4%E3%83%B3.pdf>

and

<https://www.bio.fasmac.co.jp/FasmacWebSystem/Download/%E3%82%B5%E3%83%B3%E3%83%97%E3%83%AB%E7%99%BA%E9%80%81%E6%96%B9%E6%B3%95%E3%81%AB%E3%81%A4%E3%81%84%E3%81%A6%E3%81%AE%E3%81%94%E6%A1%88%E5%86%85.pdf>