

Exo-SAP Protocol

Modified: 2010-01-27

For 1x reaction:

Reagent	Units	[Stock] (U/ μ L)	Vol (μ L)
Exo1	2	10	0.2
SAP	2.5	5	0.5
Sterile H ₂ O	---	---	1.3

Total Volume: 2.0 μ L

1) Combine 2 μ L of Exo-SAP reaction mix and 8 μ L of PCR product. (Total volume = 10 μ L)

2) Run in thermal cycler with program:

Temp ($^{\circ}$ C)	Time
37	30:00
80	15:00
4	5:00

3) Store in 4 $^{\circ}$ C.