

Plant and Animal Cell Lab

1. Make a wet mount of onion cells and water.
2. Observe under all three powers.
3. Take off cover slip and add 1-2 drops of iodine, put cover slip back
4. Observe under all three powers, and draw a picture of what you see under low or medium power, and show Mr. Howell what you see under medium power.

1. Make a wet mount of your cheek cells and water.
2. Observe under all three powers.
3. Take off cover slip and add 1-2 drops of methylene blue, put cover slip back
4. Observe under all three powers, and draw a picture of what you see under low or medium power and show Mr. Howell what you see under medium power.